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Table of contents	Model*	Page
Description of models / Delivery extent / Warranty		4
Detail view		6
Menu structure	standard, accu	8
Instructions for setup and putting into operation		10
Operation with natural gas / propane/butane gas		12/14
Flame setting/adjustment		14
Features (phoenix II eco)	eco	16
Operating modes	standard, accu	18
Operating mode “Hand“/”Foot“/”Sensor“: Setting of burn-time	standard, accu	20
Information on battery (phoenix II accu)	accu	20
Operating mode “Sensor“: Setting options of sensor feature	standard, accu	22
Operating mode “Settings”	standard, accu	24
Remarks and information in the display	standard, accu	26/28
Constructive safety measures		28
Tilting		28
Cleaning / Service		30
Troubleshooting		32
Technical Data		34
Accessories / Spare-parts		36

* all other information apply to all 3 models

!! General safety instructions !!

- **Switched on Safety Bunsen Burner should not be left unattended.**
- **Shut the gas supply when the schuett phoenix II is not in use (e.g. overnight).**
- **Operation only with provided power supply and original battery.**
- **Operation only in sufficiently ventilated rooms.**
- **Keep sufficient distance to inflammable materials and liquids.**
- **Make sure the connection to the gas-safety hose or to the gas cartridge adapter is tightly fixed.**
- **Note the required operating pressure of the corresponding gas sources: 20 mbar for natural gas, 50 mbar for propane/butane gas.**

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schuett phoenix II eco (3.354 202) with Foot-switch

Operating modes “Foot” and “Hand”

Start/stop flame by foot-switch or pressing a button by hand.

schuett phoenix II standard (3.355 202) with Sensor and Foot-switch

Colour display and language selection

Operating modes “Sensor”, “Foot” and “Hand”

Start/stop flame by activating sensor 1x or 2x, use foot-switch or press button by hand.

Burn-time displayed.

schuett phoenix II accu (3.356 202) with sensor, built-in high-power battery

(Foot-switch optional)

Colour display and language selection

Operating modes “Sensor”, “Foot” and “Hand”

Start/stop flame by activating sensor 1x or 2x, use foot-switch (optional) or press button by hand. Burn-time and battery status displayed. High-power battery integrated to the housing for up to 30 hours running time (ON_4 sec...OFF_12 sec).

Delivery extent

- schuett phoenix II Safety Bunsen Burner (eco, standard or accu)
- Foot-switch (eco, standard)
- Olive-shaped hose coupling
- Universal plug-in power supply, worldwide (eco, standard)
(set of exchange plugs – Euro, UK, US, Australia)
- Universal battery charger, worldwide (accu)
(set of exchange plugs – Euro, UK, US, Australia)
- Nozzle for natural gas (installed)
Nozzle for propane/butane gas (for exchange screwed to bottom plate)
- Operating instructions
- Warranty card with reply card

Warranty

The manufacturer grants a one-year-guarantee on the schuett phoenix II Safety Bunsen Burner. To establish your right for warranty coverage, send in the reply card accompanying the Warranty Card to the manufacturer within 14 days. Device faults and malfunctions resulting from improper use are not covered by the warranty.

Only original accessories and spare-parts are allowed for operation. Opening the burner housing without authorisation of the manufacturer voids all warranty claims.

The schuett phoenix II Safety Bunsen Burner is only suitable for intended operation as laboratory burner. The manufacturer assumes no liability for inappropriate applications. Please view the information on the enclosed warranty card.

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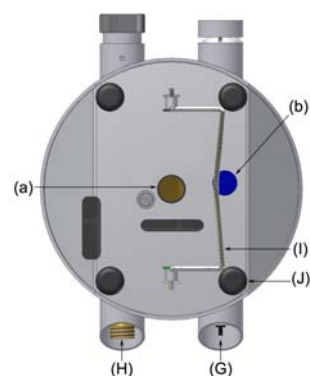
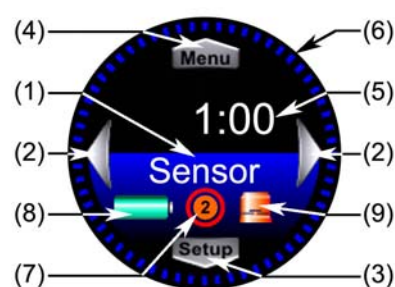
Detail view

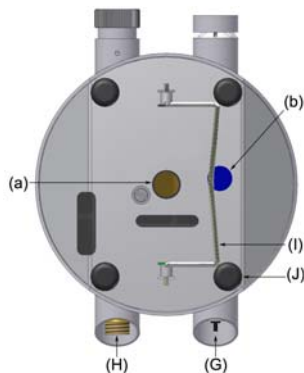
Front view from the top

- (A) Main control (Joystick)
- (B) Combined gas-air control
- (C) Sensor (standard, accu)
- (D) Display (standard, accu)
- (E) Burner head
- (F) Sparking electrode


Back view

- (G) Power connection
- (H) Gas connection (1/4" left-hand)
- (I) Bracket for tilting the burner

Bottom view

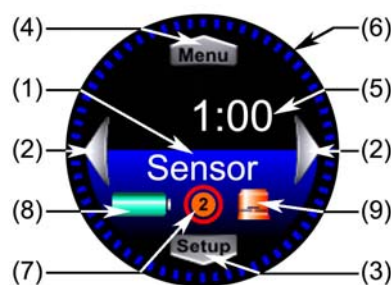- (G) Power connection
- (H) Gas connection (1/4" left-hand)
- (I) Bracket for tilting the burner
- (J) Rubber feet
- (a) Nozzle for natural gas (gold)
- (b) Nozzle for propane/butane gas (blue)


Display (standard, accu)

The display structure is consistent in all menus. The structure is shown as an example for the operating mode "Sensor" as follows:

- (1) **Display operating mode**
Hand, Foot, Sensor, Settings
- (2) **Change of operating mode**
Moving the joystick to the right or to the left
- (3) **Setup – Changing to setup levels 1-4**
Moving the joystick downwards
- (4) **Menu – Returning to setup levels or to operating modes**
Moving the joystick upwards
- (5) **Display of digital burn-time** [min:s]
Adjustable in setup level 1
- (6) **Graphical display of the digital burn-time**
Circle of seconds = 1 blue line equates to 1 s
Circle of minutes = 1 green line equates to 2 min
- (7) **Sensor activation** (only for operating mode "Sensor")
- (8) **Battery status** (only for phoenix II accu)
- (9) **Remark "Hot burner head"**

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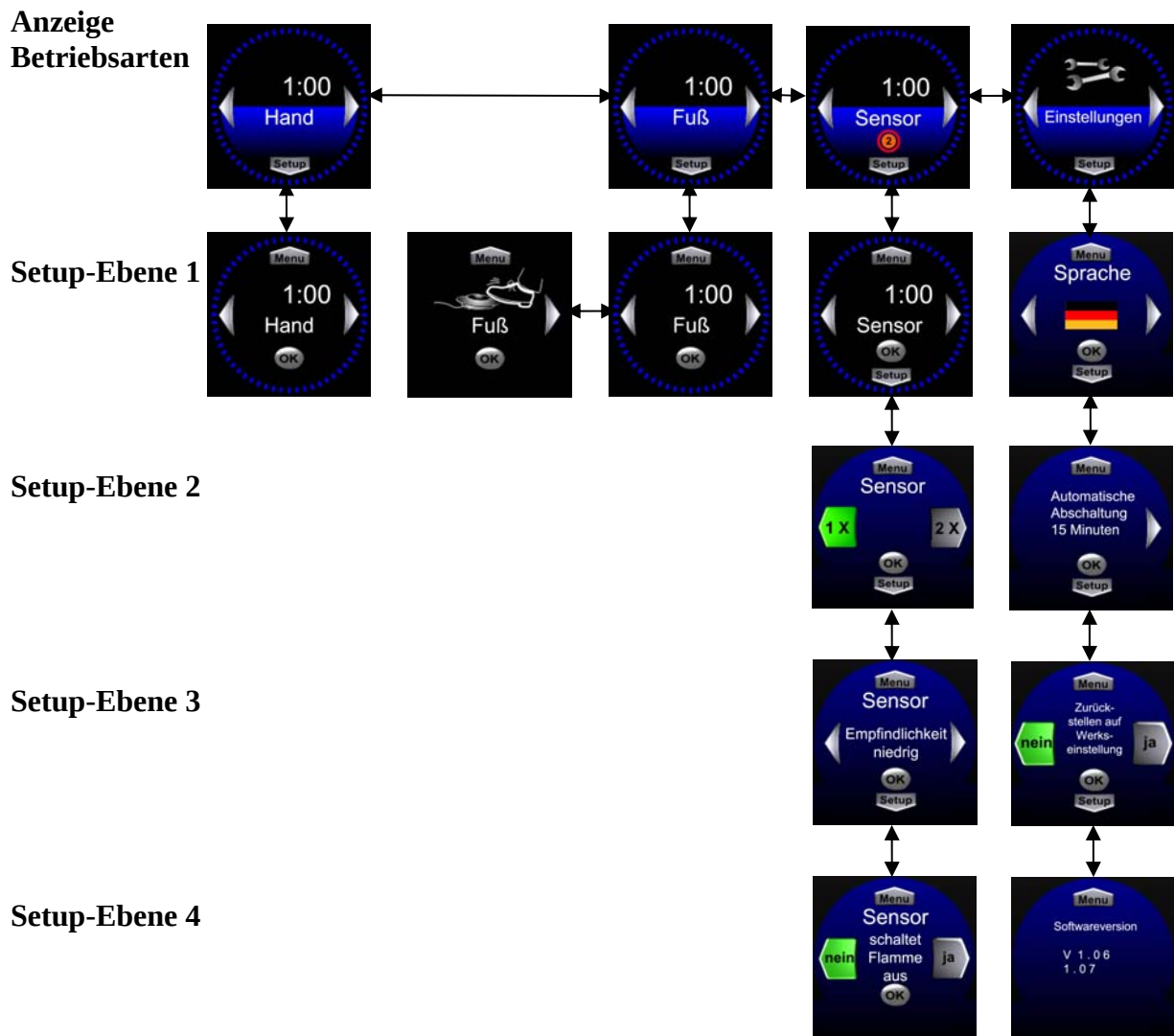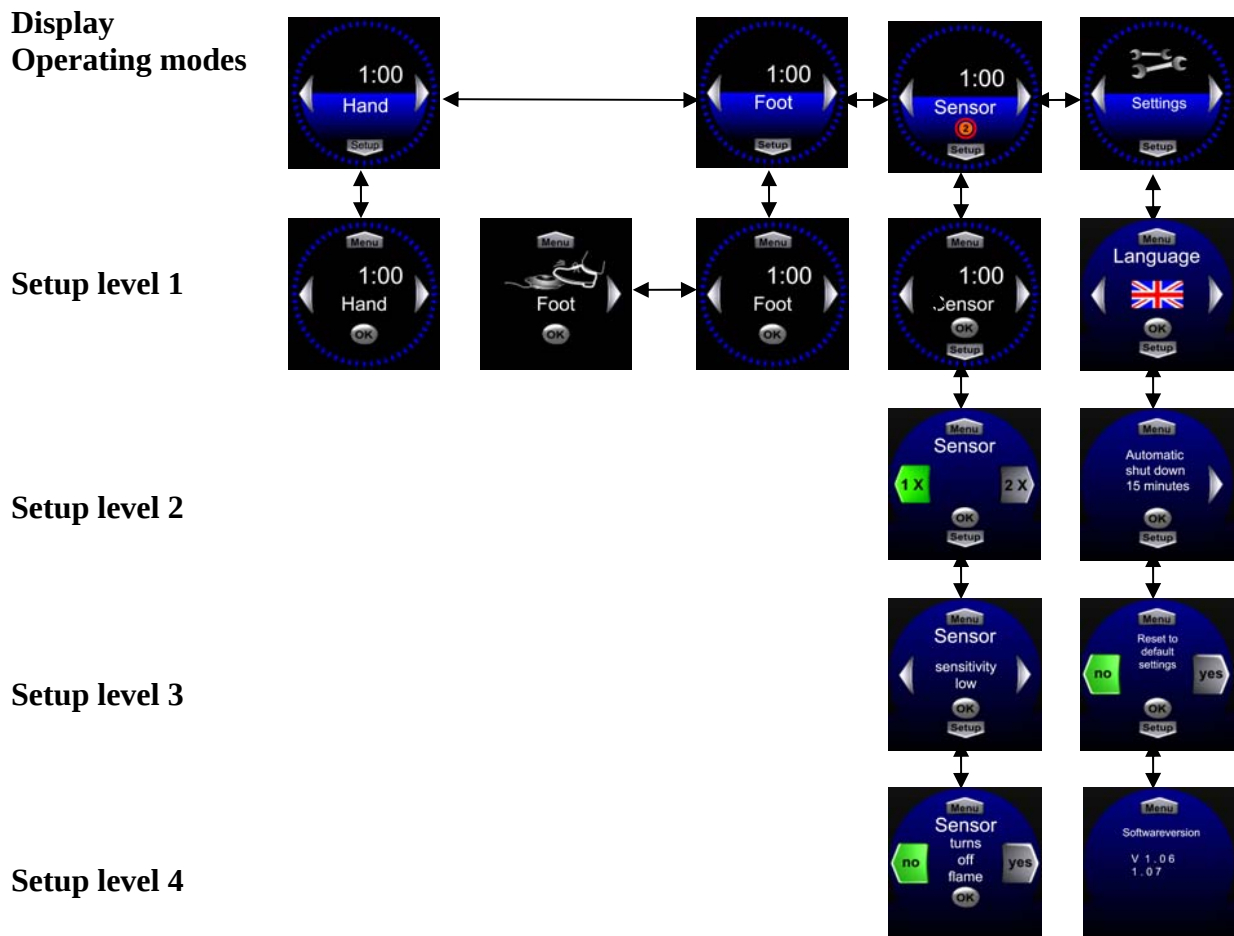
Menu structure (standard, accu)

For activating and stopping the burner, press the joystick in the centre for minimum 1 second.

↔ = Move the joystick to the right or to the left for changing the operating modes and for adjusting the settings.

↕ = Move the joystick upwards or downwards for changing between the setup levels.

 = Confirm by shortly pressing the centre of the joystick.

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Instructions for setup and putting into operation

1. Remove the device from its shipping package

Place it in its intended location on a smooth, closed surface.

Remove the transportation lock (schuett phoenix II accu only).

2. Select the gas nozzle and gas supply (for detailed information see page 12/14)

The nozzle for natural gas (gold) is mounted from the manufacturer. To operate the burner with natural gas (20 mbar operating pressure), mount the provided olive-shaped hose coupling to the **1/4"-left-hand thread** of the gas connection (H) at the burner by using a 17 mm combination wrench.

When operating the burner with propane/butane gas (50 mbar operating pressure), exchange the nozzle for natural gas by the nozzle for propane/butane gas (blue) that is screwed sideways to the centre in the housing's bottom. Mount your gas supply.

3. Connect the power supply – operation with or without foot-switch – charge battery

Compare the data on the power supply type plate with the available line (mains) voltage. The universal power supply is suitable for voltages of 110-240 V AC, 50-60 Hz, the exchange plug (Euro, UK, US, Australia) may be chosen accordingly.

a) For operation without foot-switch, connect the plug of the power supply cable with the socket (G) at the back of the burner. Please note the cut-out of the plug, the pin of the socket needs to catch the cut-out of the plug when connecting.

b) For operation with foot-switch, the power supply of the burner is **always** provided via the foot-switch. Plug the foot-switch cable to the socket (G) at the back of the burner. Connect the power supply cable to the corresponding socket on the foot-switch.

Charge the battery of the schuett phoenix II accu with the power supply for 12 hours without interruption before putting into operation for the first time (for detailed information, see page 16).

4. Control of burner head

Check whether the burner head (E) is safely caught. There is supposed to be a gap of approx. 1-2 mm between burner head and housing visible.

5. Switching the burner on

For starting the schuett phoenix II, press the centre of the joystick (A) for minimum 1 sec. The operating mode (hand) and burn-time (1:00 min) are pre-set for the first activation and are shown in the display. The burner always starts in the last chosen operating mode. A change of operating mode may be executed by moving the joystick to the right or to the left.

6. Language selection - standard, accu - (detailed information, see page 24)

The language is preset to English. Choose your language in the menu "Settings".

7. Flame setting (detailed information, see page 14)

Adjust the gas supply by moving the combined gas-air control (B) to the right or to the left and open the air supply by pushing the control fully towards the housing.

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Operation with natural gas

The nozzle for natural gas (a), golden slotted screw, is mounted in the centre of the housing's bottom upon delivery. Connect the schuett phoenix II to the gas supply by using a technically appropriate DVGW*-tested gas-safety hose (accessory). Slip the hose onto the previously mounted olive-shaped hose coupling of the gas supply (see page 10, item 2.).

Operation with propane/butane gas

The nozzle for propane/butane gas (b), blue slotted screw, is screwed sideways to the centre in the housing's bottom upon delivery. For change of nozzle, remove the nozzle for natural gas, e.g. by using a coin, and insert the required nozzle for propane/butane gas. Tighten both nozzles again. Mount the cartridge adapter or the gas-safety hose to the gas connection (H).

Operation with cartridge adapters incl. pressure reducer and gas-safety hose

Tighten the gas-safety hose's nut (**Attention: left-hand thread**) to the gas connection (H) of the schuett phoenix II by using a 17 mm combination wrench.

Make sure the stopcock of the pressure reducer is locked (turn the knurled screw to the very right).

Please note the information sheet enclosed to the adapters.

■ Cartridges without valve, e.g. C 206

Unscrew the cartridge holder from the pressure reducer and insert the cartridge. Screwing the cartridge holder onto the pressure reducer, opens the cartridge. Turn the knurled screw of the pressure reducer to the left in order to open the stopcock. The burner is now ready for operation.

Attention: When removing from the adapter, please make sure the cartridge is completely empty - gas may leak. However, the complete adapter incl. cartridge, holder and gas-safety hose may be removed from the burner at any time without gas leakage after the knurled screw of the pressure reducer has been closed at the adapter.

■ Cartridges with valve, e.g. CV 270/300/470, CG 1750, CP 250

The pressure reducer is screwed or plugged directly onto the cartridge. Open the stopcock of the pressure reducer by turning the knurled screw to the left, the burner is now ready for operation. Due to the integrated valve in the cartridge, the pressure reducer may be removed without gas leakage at any time.

Operation with CV 360-gas cartridges

Tighten the CV 360 adapter hose's nut to the gas connection thread of the schuett phoenix II (1/4"-left-hand thread) by using a 17 mm combination wrench. Make sure the stopcock of the pressure reducer is closed (knurled screw is turned to the very right). Insert the CV 360-cartridge with an audible click. In this manner, the valve in the CV 360-cartridge will be opened. Sometimes the cartridge needs to be rotated by 90° to completely open the valve. Now open the stopcock of the pressure reducer by turning the knurled screw to the left, the burner is now ready for operation. For exchanging the cartridge, simply pull the CV 360-gas cartridge out of the adapter. This locks the valve and allows for removing even partially empty cartridges from the adapter.

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Operation time in continuous mode for popular types of cartridges

When the following cartridge types are used, the schuett phoenix II achieves the following operation times for a continuously burning flame with one cartridge at 75% opened gas supply (stated operation times are only approximate values, since the operation time is also affected by regulation of the gas-flow, ambient temperature, etc.):

- CV 360-cartridge (approx. 52 g): approx. 40 minutes
- C 206-cartridge (approx. 190 g): approx. 160 minutes

Operation with 5 or 11 kg gas bottles or other cartridges

Tighten the DVGW*-certified gas-safety hose (accessory) to the gas connection thread (1/4" left-hand) of the schuett phoenix II by using a 17 mm combination wrench.

Tighten the other end of the hose to the thread of the gas supply or the pressure reducer by using a 17 mm combination wrench.

Remark: Gas bottles and cartridges have to be equipped with a DVGW*-certified pressure reducer (50 mbar) (see accessories/spare-parts, page 36).

Attention!

In case of gas odour during operation, immediately switch off the burner, close the gas supply, open a window and eliminate the cause for gas leakage.

Flame setting/adjustment

The schuett phoenix II Safety Bunsen Burner has been designed to provide a stable, sharp, hot flame when the gas and air supply are both fully opened.

Flame is too small:

Open the gas supply by turning the gas-air control (B) counter clockwise.

Flame too soft, bright:

Open the air supply by fully pushing the gas-air control towards the housing.

Flame too sharp:

In case the flame is too sharp and temporarily moves inside the burner, please decrease the air supply by pulling the gas-air control a bit out.

For a soft, bright flame:

In case of fully opened gas supply, reduce the air supply by pulling the gas-air control out. A soft, bright flame has a low temperature and is suitable for e.g. short flame-sterilizing of butyl rubber plugs or cotton plugs.

A smaller flame in total:

Decrease the gas supply. Turn the gas-air control clockwise. It is important that the air supply is reduced at the same time by pulling the gas-air control a bit out.

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Features (phoenix II eco)

Operating mode “Hand”

Activate the flame by shortly pressing the centre of the joystick.
The flame is burning until it is stopped by pressing the joystick again.

For your own safety, the maximum-burn time has been preset to 15 minutes by the manufacturer.
This setting may be increased up to 120 minutes.

At the bottom of the burner, there is a long hole under which a rotary encoder switch is situated.
Move this switch by using a small screw driver in order to adjust the desired maximum burn-time:

0 = 15 min; 1 = 30 min; 2 = 60 min; 3 = 80 min; 4 = 120 min

5 = 15 min; 6 = 30 min; 7 = 60 min; 8 = 80 min; 9 = 120 min

Operating mode “Foot-switch” (connection, see Putting into operation (page 10))

Activate the flame by shortly pressing the foot-switch. The burner is only activated as long as the foot-switch is pressed.

Indicator lamp “Power”

Red = **Burner is ON**, normal operation.

Flashing red/yellow = **Warnings** - Please check the burner.
For safety reasons the burner may not be re-activated while this remark is shown.

1. **Clean burner head:** If the flame stops due to e.g. spilled liquids and the burner may not be re-ignited, the gas supply is automatically and permanently shut. The burner head is soiled. Clean the burner head (see “cleaning the burner head”, page 30) and confirm the setting by shortly pressing the centre of the joystick.
2. **Check gas supply:** In case during automatic (re-)ignition the burner may not produce a flame after 10 seconds, the schuett phoenix II eco automatically shuts the gas supply. Please check the gas supply and confirm the setting by shortly pressing the centre of the joystick.
3. **Burner overheated:** The schuett phoenix II is suitable for continuous operation (120 min) and may not overheat. However, if the burner gets too hot due to inappropriate conditions such as back reflections of heat or unfavourable setting of flame (gas and air supply), the device automatically switches off the flame and shuts the gas supply. Wait until the burner has cooled down, the device automatically returns to the normal operating mode.

Flashing red = **Hot burner head – Caution: Risk of burns!**
After switching the schuett phoenix II eco off, the indicator lamp flashes as long as the burner head is hot. As soon as the burner head has cooled down, the flashing stops automatically.

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Operating modes (standard, accu)

Operating mode “Hand”

Activate the flame by shortly pressing the centre of the joystick.
The flame is burning until the preset burn-time has elapsed.
For interruption, shortly press the centre of the joystick again.

Operating mode “Foot” with burn-time display or foot-switch activation

Connect the foot-switch (see “Putting into operation”, page 10).

Burn-time display

Activate the flame by shortly pressing the foot-switch.
The flame is burning until the preset burn-time has elapsed.
For interruption, shortly press the foot-switch or the centre of the joystick.

Foot-switch activation

Reduce the burn-time to minimum, so that the foot-switch icon is displayed. Now, the burner is only activated as long as the foot-switch is pressed.

In case no foot-switch has been connected, the burner shows the following remark when choosing the operating mode “Foot”: “Plug in foot-switch“. Please connect the foot-switch (see “putting into operation”, page 10).

Operating mode “Sensor”

Activate the burner in sensor mode by moving quickly in front of the sensor.

For your own safety, please regard the setting opportunities for the sensor under Sensor “Setting options of sensor feature/Sensor sensitivity”, see page 22.

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Setting of burn-time (standard, accu)

Choose the requested operating mode “Hand” – “Foot” – “Sensor”.
Move the joystick downwards and choose the setup level 1.

The display shows the current preset burn-time flashing (e.g. “1:00” = 1 minute) and may be adjusted by moving the joystick to the right (+) or to the left (-).

Up to 3 minutes, the burn-time display is adjusted in seconds and after that in minutes. For rapid change of time hold the joystick to the desired side, for fine adjustment shortly move the joystick to the right or to the left.

The circle of seconds (blue) and the circle of minutes (green) visually display the burn-time in addition. One blue line equals to one second, one green line equals to two minutes.

Confirm the settings by shortly pressing the centre of the joystick,  lights up in green colour.
Return to the display of operating mode by moving the joystick upwards.

Remark:

The burner automatically switches off after the preset time, if it is no longer in use (more information under operating mode “Settings”, see page 24).

Information on battery (phoenix II accu)

- **Charge the battery of the schuett phoenix II accu with the power supply for 12 hours without interruption before putting into operation for the first time.**
- The battery charger is integrated in the power supply and provides all safety appliances. Overcharging the built-in battery is prevented.
- Charge the battery after operation or when the display shows “Charge battery”.
- Only use the built-in battery or replacement batteries from the manufacturer.
- Never use a battery that is damaged or exhausted.
- Do not expose the charging and battery contacts to conductive materials.
- Dispose exhausted batteries according to governmental regulations. Forward them to recycling. Do not throw the battery into fire.

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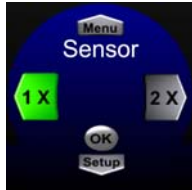
Setting options of sensor features (standard, accu)

Go to operating mode “Sensor”.

Move the joystick downwards and choose the corresponding setup level.

Setup level 2 – Setting “Sensor activation” (1x operation or 2x operation)

The activation of the sensor is adjustable.

1x operation = one movement in front of the sensor activates the flame.

2x operation = two succeeding movements in front of sensor activate the flame.

1. Move the joystick to the right, if you wish to choose “2x operation“ or to the left for “1x operation“, the arrow is highlighted in green colour.
2. Confirm the setting by shortly pressing the centre of the joystick,  lights up in green colour.

Sensor indicator lamp for 1x operation

= Burner is waiting for impulse

= Ignition and operation

Sensor indicator lamp for 2x operation

= Burner is waiting for impulse

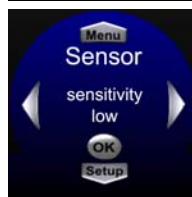
= First impulse has been recognized

= Ignition and operation

Setup level 3 – Setting of “Sensor sensitivity”

off: Sensor is switched off.

low: Sensor is activated by touching the sensor field.

medium: Sensor responds up to a distance of approx. 3 cm.

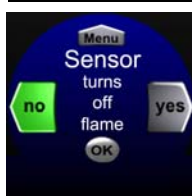
high: Sensor responds up to a distance of approx. 8 cm.
 In this position, the sensor even responds to moving an e.g. inoculating loop or glass pipette in front of the sensor field.

1. Move the joystick to the right or left until the requested sensitivity is displayed.
2. Confirm the setting by shortly pressing the centre of the joystick,  lights up in green colour.

For setting “off“, the burner shows the following remark when operating mode “Sensor” has been chosen:

Switch
on
sensor

Setup Level 4 – Setting “Sensor turns off flame”

For turning off the flame with a movement.

1. Move the joystick to the right, if you wish to turn off the flame with the sensor feature or move it to the left. The corresponding arrow is highlighted in green colour.
2. Confirm the setting by shortly pressing the centre of the joystick,  lights up in green colour.

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Setup Level 1 “Language“

1. Move the joystick to the right or left until the centre of the display shows the flag of your favoured language.
2. Confirm the setting by shortly pressing the centre of the joystick,  lights up in green colour.

Setup Level 2 “Automatic shut down”

Choose the time after which the device automatically switches off, if it is not operated.

1. You may choose between 15 min, 30 min, 60 min or 120 min.
Move the joystick to the right (+) or left (-) to set the time.
2. Confirm the setting by shortly pressing the centre of the joystick,  lights up in green colour.

Setup Level 3 “Reset to default settings”

This resets all settings to standard values (see technical data).

1. Move the joystick to the right, if you wish to reset to default settings or move to the left. The corresponding arrow is highlighted in green colour.
2. Confirm the setting by shortly pressing the centre of the joystick,  lights up in green colour.

Setup Level 4 “Softwareversion”

In case of a defect of the burner, please mention the two version numbers stated in this setup level.

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Remarks and information in the display (standard, accu)

Ignition control “Check gas supply”

In case during automatic (re-)ignition the burner may not produce a flame after 10 seconds, the schuett phoenix II automatically and permanently shuts the gas supply. The display shows “Check gas supply“. See “Troubleshooting”, page 32.

Permanent burning control “Clean burner head”

If the flame stops due to e.g. spilled liquids and the burner may not be re-ignited, the gas supply is automatically and permanently shut. The burner head is soiled and the display shows the remark “Clean burner head”.
See “Troubleshooting”, page 32.

Warning “Hot burner head”

If the flame has been burning for a longer time, the burner head is hot, the display mentions this . Even after switching the burner off, a blinking remark is shown in the display. It disappears automatically as soon as the burner head has cooled down.

Wait until this remark vanishes, touching the burner head before may cause burns!

Warning “Burner overheated”

The schuett phoenix II is suitable for continuous operation (120 min) at fully opened gas and air supply, the burner may not overheat. However, if the burner gets too hot due to inappropriate conditions such as back reflections of heat or unfavourable setting of flame, the device automatically switches off. The remark “Burner overheated” is shown in the display.

Wait until the burner has cooled down, the display automatically returns to the last chosen operating mode.

Controlled burn-time/remaining time – safety shutdown

Digital and exact to the second adjustable (1 sec ... 120 min), safety shutdown after 120 minutes.
See “Settings automatic shutdown”, page 24.

Sensor distance (approx. 0 to 80 mm)

The distance is adjustable in 4 steps.

See “Setting options of sensor features/Setting of sensor sensitivity”, page 22.

Adjustable sensor activation

1x operation /2x operation, display of sensor indicator lamps, sensor turns flame off.

See “Setting options of sensor features/Setting of sensor sensitivity”, page 22.

Flame control

In case of unexpected extinguishing of the flame, it is automatically re-ignited, so that the work does not need to be interrupted.

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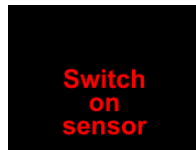
“Switch on sensor”

In case the sensor sensitivity is set to “off”, the burner shows the following remark shortly after the operating mode “Sensor” has been selected.
See “Setting options of sensor features/Setting of sensor sensitivity”, page 22.

„Plug in foot-switch“

If the foot-switch is not connected and the operating mode “Foot” is selected, the burner shows the remark “Plug in foot-switch” shortly afterwards.
Please connect the foot-switch (see “Putting into operation”, page 10).

“Charge battery” (only for battery operation)

The battery has fallen below its minimum voltage. Charge the battery by using the provided battery charger at the line/mains supply.

Certified safety:

The device meets the latest safety regulations such as DIN 30665 chapter 1 and has been registered under DIN-DVGW* certificate NG-2211AR0750.

Constructive safety measures

Low placed burner tube and sparking electrode

The burner tube and sparking electrode are placed low inside the burner head to guarantee a stable flame, even in strong currents in laminar flow cabinets.

Drain for liquids

Liquids entering the burner head are drained via a special opening in the bottom. Clogging of the air regulation is not possible due to its advanced design.

Gas Input filter

A filter in the magnet valve prevents particles from entering the burner via the gas supply.

Overheating protection

The burner head of the schuett phoenix II is physically separated from the burner by a thermally insulating jacket, i.e. even in the continuous mode the temperature of the housing never exceeds 60 °C.

Tilting

For flame-sterilizing e.g. bottlenecks, it can be advantageous to put the schuett phoenix II in a tilted position. Switch the burner off and lift the bracket until it locks in place. Tilt the burner accordingly, depending on left-hand or right-hand operation.

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Cleaning

Cleaning the Safety Bunsen Burner

Clean the schuett phoenix II only with commercial cleaning solutions and soft cloths. The smooth stainless steel and tempered glass surfaces are resistant to ethanol and disinfectants. The stainless steel cover may shortly be flame-sterilized without any problems.

Remove visible soiling upon each operation and keep the display as well as the two sensor fields clean.

Cleaning the burner head

The burner head may be removed for cleaning or replacement in case of defects (see accessories/spare-parts, page 36).

Please clean the burner head if liquids were accidentally spilled over the burner. Remove the burner head and clean it in a standard dishwasher. Dry the burner head in a drying cabinet afterwards.

The burner head is attached to the burner with a bayonet catch and may therefore be demounted without tools.

- For demounting the burner head carefully turn it clockwise as far as it will go and pull it straight up out of the burner.
- For mounting, insert the burner tube of the burner head into the insulating piece of the burner, which is only possible in one position, and carefully turn the burner head counter clockwise as far as it will go. The sparking electrode is facing the back of the burner. The gap between the correctly mounted burner head and the housing is 1-2 mm.

Service

The schuett phoenix II Safety Bunsen Burner is generally maintenance-free. Please check the gas-safety hose regularly for damages and replace it, if need is.

Furthermore we advise an annual inspection of the burner:

- Check the seal of the nozzles for cracks.
- Check the burner head for soiling and make sure, the sparking electrode is fixed. Remove and mount the burner head or replace it, if need is.
- Check all cables (foot-switch/ power cable) and gas-safety hoses for damages.

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Troubleshooting

A simple fault may often be the reason for a malfunction, and as such can easily be corrected by the user himself.

Malfunction	Measures to eliminate the fault
Blank display after pressing the centre of the joystick.	▪ Check line/mains connections and voltage. ▪ Check cable connections. ▪ Is the power supply provided via foot-switch, loosen the plug and connect the line/mains cable directly to the schuett phoenix II.
After switching the burner on, it tries to ignite (audible clicking) for approx. 10 seconds, but no flame is produced. The display shows the remark "Check gas supply".	▪ Fully open the gas supply "Gas" (B) by turning counter clockwise. ▪ Check all intermediary gas taps. ▪ Replace empty cartridge or bottle. ▪ Press the centre of the joystick (A) in the hand operating mode and start the burner again.
The flame is too large when using propane/butane.	▪ Exchange the nozzle for natural gas (gold) with the nozzle for propane/butane (blue).
The flame is too small when using natural gas.	▪ Exchange the nozzle for propane/butane (blue) with the nozzle for natural gas (gold).
The flame is small or unstable.	▪ Optimize the relation of gas/air mixture by turning and pulling the gas-air control (B).
The IR-sensor does not work.	▪ Check the sensor settings (see page 22). ▪ Activate the sensor. ▪ Increase the sensitivity of the sensor.
The foot-switch does not work.	▪ Check cable connections. ▪ Check the 5 pins in the plug. ▪ If the burner works in sensor mode, exchange the foot-switch (accessory).
After switching the burner on, no ignition (no audible clicking), no flame produced. After approx. 10 seconds, the remark "Clean burner head" is shown in the display.	▪ Clean the burner head as described (see page 30) and mount the clean burner head to the device. ▪ Press the centre of the joystick (A) shortly and activate the burner again. If this does not correct the fault, please replace the burner head (spare-part).

If the measures listed above do not restore the device to its operational state or if the fault is not listed above, please contact the manufacturer.

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Technical data

Burn-time, adjustable (standard/accu)	1 sec...120 min
Line/mains voltage for the power supply	110 - 240 V AC, +/- 5%, 50-60 Hz
Connected voltage (standard/accu)	6 V / DC, 3 VA max.
Olive-shaped hose coupling suitable for natural gas hose	10 mm diameter (inner)
Working pressure for natural gas	20 mbar
Working pressure for propane/butane	50 mbar
Rated heat load, natural gas	600 W
Rated heat load, propane/butane	1,200 W
Consumption natural gas	0.06 cbm/h
Consumption propane/butane	77.2 g/h
Dimensions (W x H x D)	128 x 85 x 160 mm (incl. burner head)
Weight	approx. 1,000 g

Default settings

Language	English
Operating mode	Hand
Automatic shut down of the burner	after 120 min

Burn-time:

Operating mode "Hand"	5 min
Operating mode "Foot"	1 min 30 sec
Operating mode "Sensor"	3 sec

Sensor settings

Sensor activation	1x operation
Sensor sensitivity	Low
Sensor turns off flame	No

Operating time in battery mode (only phoenix II accu):

Continuous mode	approx. 15 h
Flame-sterilization of inoculating loops (ON_4 sec...OFF_12 sec)	approx. 30 h

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Accessories (optional) / Spare-parts	Cat.-no.
Foot-switch (works in 2 operation modes) 3500 mm cable length**	3.358 392
Glass spatter guard, to work with pathogene materials	3.358 332
Windshield stainless steel, to protect the flame from sideways currents	3.358 342
Instrument Tray stainless steel, with 5 holding spaces	3.358 352
Hot-Tray stainless steel, for use in chemical laboratories, height-optimized	3.358 362
Adapter cable for connection of 12 V (car cigarette lighter)	3.358 382
Pressure reducer for 5 kg and 11 kg propane gas bottles (TÜV-certified, indoor/outdoor)	3.358 302
Gas-safety hose for natural gas with standard plug 1/2" (1,0 m / 2,0 m)** for the connection to a gas-outlet with thermal shut-off device (TAE)	3.358 512 / 3.358 522
Burner head (standard)	3.935 700
Burner head (long)	3.935 710
Universal plug-in power supply (eco/standard), 100-240 V, 50/60 Hz, worldwide applicable (incl. set of exchange plugs – Euro, UK, US, Australia)	3.935 800
Universal battery charger (accu), 100-240 V, 50/60 Hz, worldwide applicable (incl. set of exchange plugs – Euro, UK, US, Australia)	3.935 810
Spare-part nozzle for propane/butane gas	3.935 730
Spare-part nozzle for natural gas	3.935 740
Spare-part olive-shaped hose coupling	3.935 790
Adapter for natural gas connection 1/2" (France)	3.934 192
Gas-safety hoses	
for natural gas (0.75 m / 1.5 m / 2.0 m)	3.359 212 / 3.359 222 / 3.359 232
for propane/butane gas, with threaded connections (0.5 m / 2.0 m)	3.359 272 / 3.359 282
Adapter incl. pressure reducer for connection	
of gas cartridges CV 360, incl. pressure reducer	3.357 212
of gas cartridges Express 444	3.357 362
Adapter incl. pressure reducer and gas-safety hose (0.5m) for connection	
of gas cartridges C 206	3.357 232
of gas cartridges CV 270 / CV 300 / CV 470	3.357 252
of gas cartridges CP 250	3.357 302
of gas cartridges CG 1750	3.357 322
of gas cartridges CG 901 / CG 904 / CG 907	3.357 342
Gas cartridges	
Cartridge CV 360, with valve, 52 g/cartridge butane gas (qty=6)	3.351 300
Cartridge C 206, without valve (pierceable), 190 g/cartridge butane gas (qty =6)	3.351 310
Cartridge CV 470, with valve, 450 g/cartridge butane gas (qty =3)	3.351 342
Cartridge CG 1750, with screw-valve, 175 g/cartridge propane/butane gas (qty =6)	3.351 352
Cartridge CP 250, with valve, 250 g/cartridge butane gas (qty =6)	3.351 362
Cartridge CV 270, with valve, 230 g/cartridge propane/butane gas (qty =6)	3.351 372
Cartridge CV 300, with valve, 240 g/cartridge propane/butane gas (qty =6)	3.351 392
Cartridge Express 444, with valve, 60 g/cartridge propane/butane gas (qty=6) (GGVSEB 2/5 F, UN 2037, LQ 2)	3.351 382

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schuett-biotec GmbH